



Code 2984LR


Peroxide Low Range
25 Test Strips



Insta-Test[®] ANALYTIC

Peroxide ppm	Code 2984LR
0	
1	
3	
10	
30	
50	

Using a cup-size sample, immerse strip for **2 seconds**. Remove strip with **pad face up**. **DO NOT SHAKE OFF EXCESS WATER**. With pad still facing up, immediately compare to color chart. Discard strip after 30 seconds.

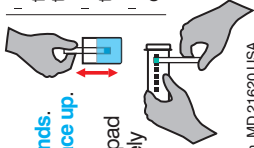

- Keep wet fingers out of the bottle.
- Cap bottle tightly.
- Store in a cool, dry place.



LaMotte Company
802 Washington Ave., Chestertown, MD 21620 USA
410-778-3100 • www.lamotte.com

