

TOTAL HARDNESS KIT

DROP COUNT, 1 DROP = 2, 5, OR 10 ppm

CODE 7246-02

QUANTITY	CONTENTS	CODE
2 x 30 mL	*Hardness Reagent #5	*4483-G
30 mL	*Hardness Reagent #6 Solution	*4485-G
60 mL	Hardness Titrant 2	2788WT-H
1	Sample Tube, 5-10-25 mL, w/cap	0715

*WARNING: Reagents marked with an * are considered to be potential health hazards. To view or print a Safety Data Sheet (SDS) for these reagents go to www.lamotte.com. Search for the four digit reagent code number listed on the reagent label, in the contents list or in the test procedures. Omit any letter that follows or precedes the four digit code number. For example, if the code is 4450WT-H, search 4450. To obtain a printed copy, contact LaMotte by email, phone or fax.

Emergency information for all LaMotte reagents is available from Chem-Tel: (US, 1-800-255-3924) (International, call collect, 813-248-0585).

To order individual reagents or test kit components, use the specified code number.

NOTE: This test allows the analyst to use different sample volumes to vary equivalencies. Select the appropriate sample volume from the table below and add reagents as specified.

SAMPLE SIZE	EQUIVALENCE (ppm per drop)
25 mL	1 drop = 2 ppm
10 mL	1 drop = 5 ppm
5 mL	1 drop = 10 ppm



Yalitech Instruments

Santiago de Chile, Río Refugio 9648, Parque de Negocios ENEA, Pudahuel.

www.yalitech.cl · ventas@yalitech.cl · (+56 2) 28988221

PROCEDURE

1. Rinse sample tube with water to be tested. Fill with desired sample size.
2. Add *Hardness Reagent #5 (4483) as follows:

25 mL sample	10 drops
10 mL sample	8 drops
5 mL sample	5 drops

Swirl to mix.

3. Add *Hardness Reagent #6 Solution (4485) as follows:

25 mL sample	5 drops
10 mL sample	3 drops
5 mL sample	2 drops

Swirl to mix. Solution will turn red if hardness is present. If solution is blue, there is no measurable amount of hardness.

4. Add Hardness Titrant 2 (2788WT), one drop at a time while swirling and counting the drops, until red color turns to blue. Hold bottle vertically.
5. Multiply number of drops used in Step 4 as follows:

25 mL sample	multiply by 2
10 mL sample	multiply by 5
5 mL sample	multiply by 10

Record as ppm Hardness as CaCO_3 .

LaMOTTE COMPANY

Helping People Solve Analytical Challenges

PO Box 329 · Chestertown · Maryland · 21620 · USA
800-344-3100 · 410-778-3100 [Outside U.S.A.] · Fax 410-778-6394
Visit us on the web at www.lamotte.com